

Applications

- Photovoltaic Testing
- Environmental Testing
- Photobiology
- Photochemistry
- Material and degradation testing

Features

- Class AAA specification (ASTM, IEC)
- Illumination area: 50x50mm
- Touchscreen power supply with control software included
- Manual shutter included (electronic shutter available)
- Variable attenuator from 0.1–2 suns
- Plug and play operation
- Long working distance can facilitate glovebox integration

Small Area Solar Simulators SciSun Series



SciSun Solar Simulators

OVERVIEW

Sciencetech's line of SciSun solar simulators are easy to use, economically priced, and technically superior. The SciSun line is designed for researchers who do not require a large field of illumination. They can produce up to **2 Suns** and feature Class AAA specifications.

The SciSun series provides a flexible output orientation that can be adapted to different requirements. The standard configuration is downward-facing; however, a horizontal output can be achieved easily.

Our **SciSun-LP** is an affordable Model, which include a basic power supply, and does not include some features of our Standard Model (See right table).

| Configuration | Standard Model | LP Model |
|---|----------------|----------|
| Arc Lamp Housing with integrated igniter | ● | ● |
| Xenon arc lamp | ● | ● |
| Filter holder | ● | ● |
| Beam turner | ● | ● |
| Continuous beam angle variations from 0-360 degrees | ● | ● |
| Quality Control Report | ● | ● |
| Basic Power Supply | | ● |
| Touchscreen power supply interface | ● | |
| Power supply control software | ● | |
| Manual variable attenuator | ● | |
| Height adjustable stand | ● | |

| Model | SciSun-300 | SciSun-LP-300 | SciSun-150 | SciSun-LP-150 |
|--|---|---------------|----------------------------|---------------|
| | 160-9101 | 160-9104 | 160-9103 | 160-9105 |
| Target Area | 50 × 50 mm | | | |
| Solar Simulator Class | A-A-A (Irradiance Uniformity ¹ , Spectral Match ² , Temporal Stability ³) | | | |
| Irradiance at Target (AM1.5G 1 Sun=100mW/cm ²) | Up to 2 Sun ² | | Up to 1 Sun ² | |
| Lamp Wattage (watts) | 300 | | 150 | |
| Lamp Type | Xenon Short Arc , Ozone free | | | |
| Working Distance (mm) | 380 ± 15 | | | |
| Manual Shutter | Included | | | |
| Manual Variable Attenuator | Included | Available | Included | Available |
| Dimensions (L×W×H) | 535 × 183 × 188 mm | | | |
| Weight without PS (kg) | 8.5 + 8 (stand) | 8.5 | 8.5 + 8 (stand) | 8.5 |
| Power Supply Model | 601-300 | EPS-300 | 601-150 | EPS-150 |
| Power Requirements | 110-240V, 50Hz/60Hz , 450W | | 110-240V, 50Hz/60Hz , 250W | |
| Stability / Ripple / Regulation | 0.05% / < 1% / 0.02% current variation for 5V line charge | | | |

STANDARDS

SciSun solar simulator specifications listed are according to
 ASTM E927-19
 IEC-60904-9
 JIS C8904-9 (2017)



1) Determined from true Isc measurements with silicon sensor mounted on 2 axis automated stage. 2) Measured using NIST traceable secondary reference cell. 3) Measured with scanning spectroradiometer calibrated as per ASTM G138-06. 4) Determined from 20 measurements spaced at 250ms, NPLC=1. Due to our continuous improvement system, all specifications are subject to change without notice.

SciSun Solar Simulators

CLASSIFICATION AAA

Class A Spectral Match

SciSun solar simulators match Class A spectral match when used with a compatible air mass filter (sold separately; see below using an AM1.5G filter). All testing results are for an example SciSun-300 and individual reports will vary

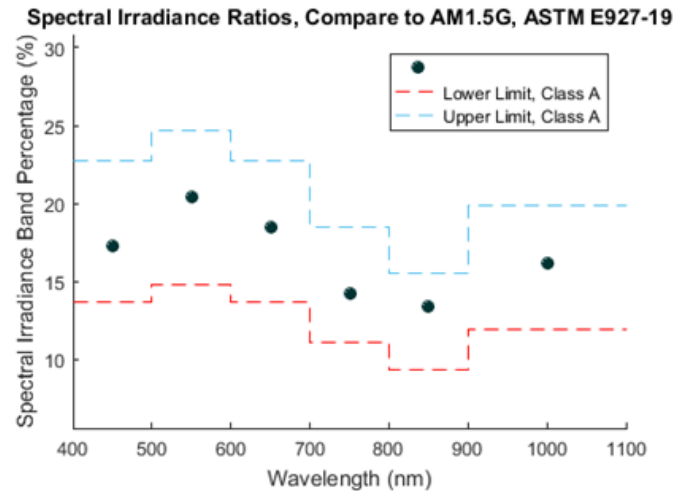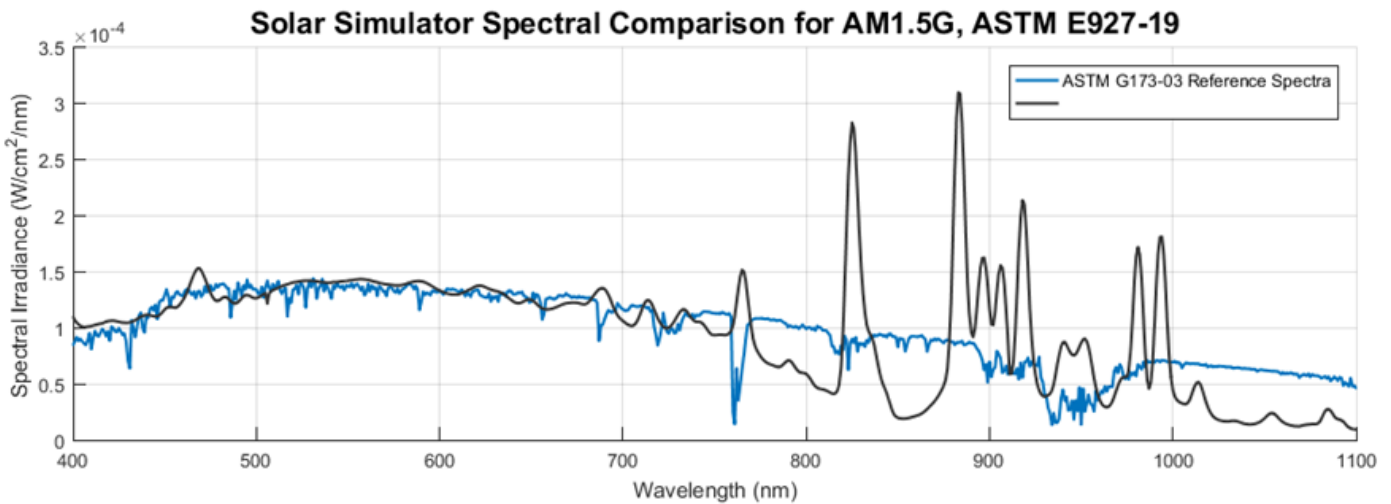

| Wavelength | Percentage | Class |
|------------|------------|-------|
| '400-500' | '17.2821' | 'A' |
| '500-600' | '20.4084' | 'A' |
| '600-700' | '18.5142' | 'A' |
| '700-800' | '14.2414' | 'A' |
| '800-900' | '13.4281' | 'A' |
| '900-1100' | '16.1258' | 'A' |

Solar Simulator Standards

SciSun solar simulator specifications listed are according to ASTM E927-19 and IEC-60904-9 unless otherwise stated. We can accommodate testing to match several standards. Testing procedure as per ASTM E927-19 provided by default. Please specify upon ordering if testing against IEC-60904-9 is required.

SciSun Solar Simulators


CLASSIFICATION AAA

Class A spatial non-uniformity (NU):

SciSun solar simulators meet Class A spatial non-uniformity by default (see below).

Non-uniformity = 1.6% Less than 2%.

Class B may also be available over larger target sizes upon request.

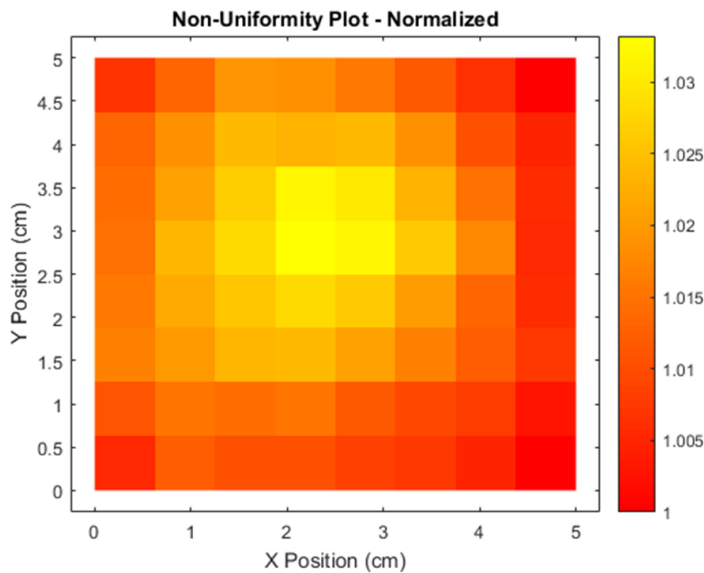


Non Uniformity Map of a SciSun Solar Simulator.

| | |
|---|----------------------|
| Detector Area: | 0.28 cm ² |
| Number of Measurement Points: | 64 |
| Measurement Point Area: | 0.39 cm ² |
| Maximum Irradiance: | 1.0169 Suns |
| Minimum Irradiance: | 0.9842 Suns |
| Sample Standard Deviation of Spatial Non-Uniformity: | 0.008 Suns |
| Spatial Non-Uniformity of Irradiance: | 1.60% |
| Classification: | A |

Class A Temporal Instability:

SciSun solar simulators meet Class A temporal instability. 0.05% Less than 2%.

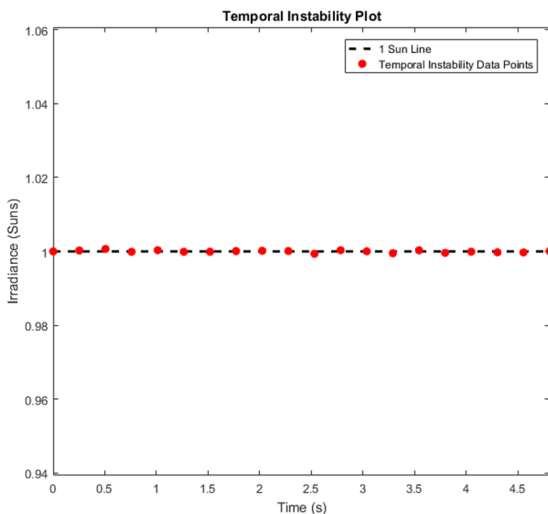

| | |
|--|-------------------|
| Detector Area: | 4 cm ² |
| Time Between Data Points: | 0.253 Sec |
| Number of Power Line Cycles (NPLC): | 1 |
| Total Measurement Points: | 20 |
| Maximum Irradiance: | 1.0007 Suns |
| Minimum Irradiance: | 0.9994 Suns |
| Temporal Instability of Irradiance: | 0.05% |

SciSun Solar Simulators

STANDARD FEATURES

Touchscreen Power Supply - 601

Each SciSun series solar simulator (non-LP series) comes with a 601-series power supply.



601-series power supply

Standard features included with Sciencetech's 601-series power supplies:

- Touchscreen interface
- Shutter and exposure control (if electronic shutter is supplied)
- Single connection for lamp power, cooling, and communication
- Lamp starts and timer log
- Fan cooling safety interlock
- RS232 software GUI included

Filter Box Assembly


This system has a modular optics assembly which can hold a range of filters in Sciencetech's standard FT style filter holder. The most popular options are AM filters; however a range of other filter options are available such as bandpass filters and neutral density filters.

Variable Aperture


Sciencetech's SciSun solar simulators include a variable aperture component, which allows variation of the output irradiance level without adjusting the power supply. The range of attenuation is continuously variable from 10% to 100%. Uniformity is best maintained at specific output levels. Non-uniformity versus output level for the VAR-ATTN-M may vary between models.

The SciSun-LP series features the EPS-series simplified power supply, which lacks a touchscreen, computer control, or control of electronic accessories.

601-series touch screen power supply main control screen

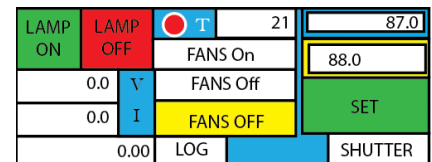


601-series touchscreen power supply automatic shutter control screen

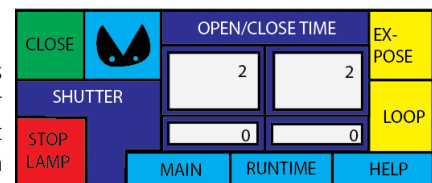

Software Included

SciSun Solar Simulator come with SciLampPower Control.

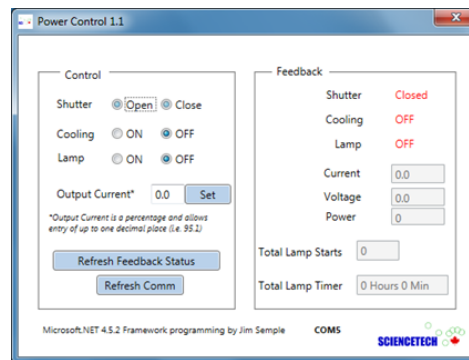

Software GUI for power supply control

Spectral Filter Options

AM1.5-FT-3 Model: Includes AM1.5 Filter-Class A with 160-8085 part number.

AM1.0D-FT-3 Model: Includes AM1.0D Filter-Class A with 160-8086 part number.

AM1.5D-FT-3 Model: Includes AM1.5 Filter-Class A with 160-8087 part number.

SciSun Solar Simulators

ACCESSORIES

**SOL-METER****(125-9011)**

Solar Power Meter, a digital meter for use with solar calibrated detectors (e.g. SSIVT-REF or SC-LT-Q).

**SSIVT-REF****(125-9007)**

This silicon detector is designed to be used for monitoring and verifying the sun level of solar simulators.

**SC-LT-Q****(585-0154)**

Calibrated Reference Cell, Quartz Window, traceable to NIST and NREL.

**SSIVT-20C****(175-9103)**

20W IV Tester for Continuous Solar Simulators (current range = 1 A, voltage range = 200 V).

**SCI-SCC3-TE****(165-8202)**

3.5" x 3.5" Solar Cell Chuck, TE Cooled, Computer controllable, Vacuum Ready.

**SCI-SCC3-L-B****(165-8221)**


3.5" x 3.5" Solar Cell Chuck, Liquid Cooled, Rear Contact.

SciSun Solar Simulators

ACCESSORIES

**SCP-4T****(165-8211)**

Probe Station, 4 Probes, Tungsten Needle-tip Kelvin Probes

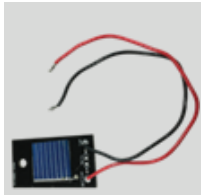**SCI-REF-NL****(125-9028)**

A simple PCB mounted solar cell, as a reference cell. No load.

**Height Adjustable Stand****(101-8024)**

Height adjustable Stand. The stand allows to adjust the working distance as required.

**MF-49-FT-3****(640-9006)**

Standard 75 mm (3") neutral density mesh filters (different light transmissions available).


**SH-SC3****(127-8004)**

Computer controlled shutter. Installs inside SciSun (non-LP versions only).


SciSun Solar Simulators

DIMENSIONS


Dimensions are in [mm] and inches.

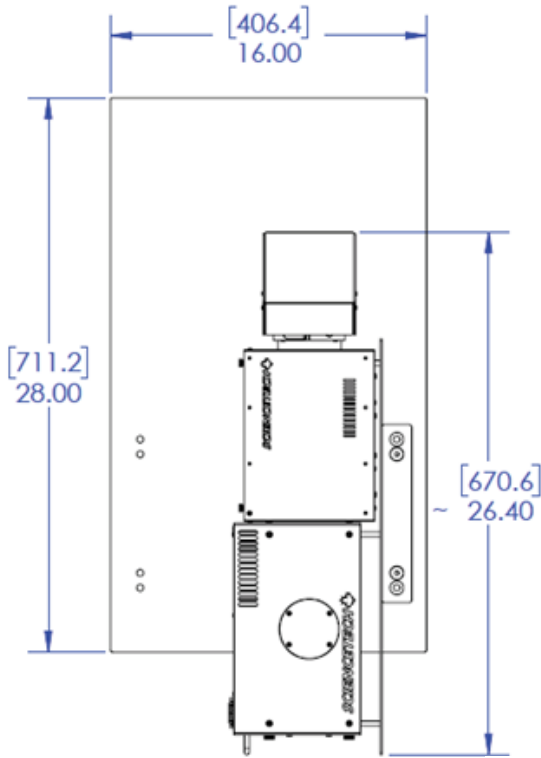

SciSun-150 and SciSun-300

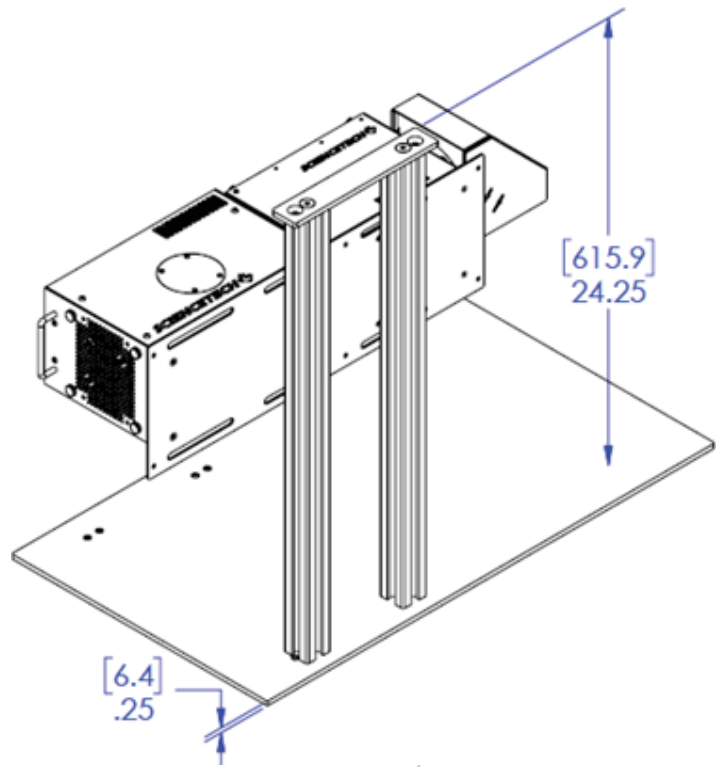

SciSun-150 and SciSun-300

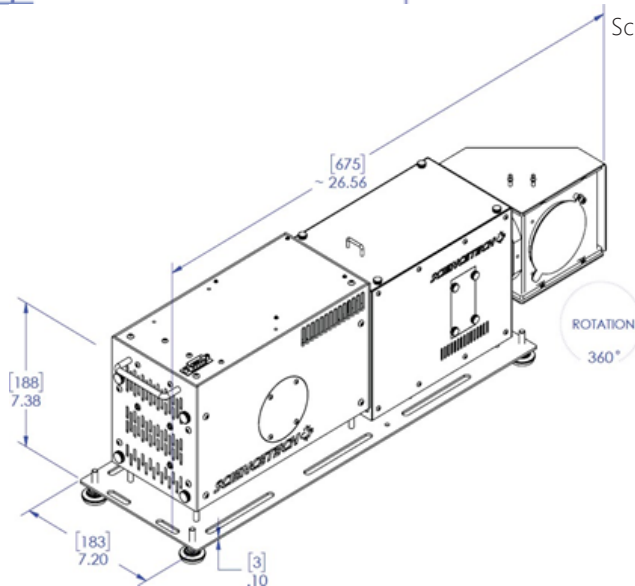

SciSun-150 and SciSun-300

Please note: Due to our continuous improvement system, all specifications are subject to change without notice. SciSun solar simulator specifications listed are according to ASTM E927 -19 and IEC-60904-9 unless otherwise stated.